(limited countries and region)





VIP ECO

Natural Refrigerants -86°C Ultra-Low Temperature Freezer









528 L / 729 L

Cost-saving and environmentally friendly sample storage within an optimal footprint.

VIP ECO -86°C Upright Freezer provides maximum sample storage capacity within an optimum footprint. When combined with natural refrigerants to minimise energy consumption, environmental impact is reduced.

Natural Refrigerants and Inverter Technology

Naturally occuring hydrocarbon [HC] refrigerants have minimal effect on the environment and are compliant with environmental legislation for climate control. Combined with inverter technology, these refrigerants also provide more efficient cooling without compromising cooling performance and ambient tolerance and recovery speeds following door openings.

Efficient Sample Storage

New cabinet design combines an insulated outer door with superior storage efficiency and an inner door designed for uniform temperature distribution. This improves storage efficiency approximately 10 % (576 pcs of 2-inch Cryo boxes can be stored in the MDF-DU702VH, MDF-DU702VHL) while maintaining the same footprint as conventional models.

Intelligent Interface

The newly developed ergonomic "EZlatch" door handle makes access to stored samples even easier. Intuitive and intelligible large colour LCD touchpanel is provided.



Energy efficient performance



Natural refrigerants, compressors and integrated electronics combine to lower operating costs by up to 30%*. Freezer operation is managed by effectively balancing temperature performance and energy management.

* Compared to conventional models. Power source: 230V/240V 50Hz, AT 30°C

Flexible shelf lavout



Reliable controllability and data log function




Large colour LCD touchpanel is accurately controlled even with a gloved hand, while the USB port makes transferring logged data of product's operational status to a PC convenient.

VIP ECO Natural Refrigerants -86°C Ultra-Low Temperature Freezer

Approximately 110 mm*1 installation width can be reduced when installing 3 units in line compared to conventional units.

The "Chamfer-design" body front features a chamfered edge for reduced installation space when compared to conventional models in a multiple unit installation. The design is especially suitable for a biobank or similar application.

*1 Comparison of MDF-DU702VH MDF-DU702VHL and the conve tional MDF-U700VH product. Comparison condition: Minimum width required when installing 3 units in line. This prevents outer doors from contacting each other for door openings/closings.

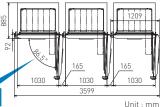


User friendly for daily operation with newly developed vacuum release port and fivefold gasket

Newly developed automatic vacuum release port and fivefold gasket are adopted. Automatic release of negative pressure and reduced frost realise user friendliness even for frequent door opening.




Model Number⁹

Performance Data

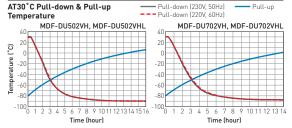

Dimensions

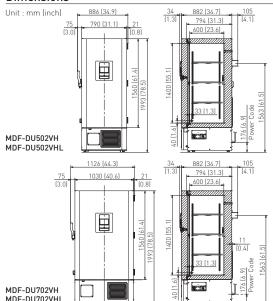

Model Number		MDF-DU502VHL-PS MDF-DU502VH-PE	MDF-DU702VHL-PS MDF-DU702VH-
External dimensions (W x D x H) 1	mm	790 x 882 x 1993	1030 x 882 x 1993
Internal dimensions (W x D x H)	mm	630 x 600 x 1400	870 x 600 x 1400
Volume	litres	528	729
Net weight	kg	246	278
Capacity 2)	2" boxes	384	576
Performance	Z BOXES	304	370
	°C		2/
Cooling performance 3		-86	
Temperature setting range	°C	-50 to -90	
Temperature control range 31	°C	-50 t	o -86
Control			
Controller		Microprocessor, non-volatile memory	
Display		LCD Touchpanel	
Temperature sensor		Pt-1000	
Refrigeration			
Refrigeration system		Cas	cade
Compressors	W	2 x 750	
Refrigerant		HC	
Insulation material		PUF / VIP PLUS	
Insulation thickness	mm	80	
Construction			
		5 : .	1.61
Exterior material		Painted Steel	
Interior material		Painted Steel	
Outer door	qty	1	
Outer door lock		Y	
Inner doors	qty	2 pieces	
Shelves	qty	3 (stainless steel)	
Max. load - per shelf	kg	50	
Max. load - total	kg	415 4) 515 4)	
Vacuum release port		2 (1 automatic, 1 manual)	
Access port	qty	3	
Access port position	17	3 locations [rear x 1, bottom x 2]	
Access port diameter	Ømm	17	
Casters		4 (2 leveling feet)	
Alarms	qty		
Power failure		(V = Visual Alarm, B = Buzzer Alarm, R = Remote Alarm)	
		V-B-R	
High temperature		V-B-R	
Low temperature		V-B-R	
Filter		V	
Door open		V-	-В
Electrical and Noise Level			
Power supply	V	PA: 220 / PE: 220/230/240 / PS: 230	
Frequency	Hz	PA: 60 / PE: 50 / PS: 50	
Noise level 5]	dB (A)	52	
Options			
	set of 5	MDF-5ID5-PW 6]	MDF-7ID5-PW 61
Small inner door kit	set of 4	MDF-5ID4-PW	MDF-7ID4-PW
Liquid CO ₂ back-up	300 01 4		
Inventory rack		MDF-UB7-PW IR-224U-PW, IR-316U-PW	
*			
e - Circular type		MTR-G85C-PE ⁷¹ - Chart paper: RP-G85-PW - Ink pen: PG-R-PW	
รับ อ Continuous strip type	·	MTR-85H-PW ⁷⁾ - Chart paper: RP-85-PW	
၅ ဗို - Continuous strip type		- Ink pen: DF-38FP-PW	
Ter			er housing: MDF-S3085-PW
Optional Communication Systems			
Digital interface (RS232C/RS485) 81		MTR-480-PW	
Ethernet interface (LAN) ⁸⁾		MTR-L03-PW	
Quality Management System			
<u> </u>		100	9001
Certification			

- 1) Exterior dimensions of main cabinet only, excluding

- 11 Exterior dimensions of main cabinet only, excluding handle and other external projections.
 21 Usable storage capacity of 2" boxes when the original 3 trays installed.
 31 Air temperature measured at freezer centre, ambient temperature 30°C, no load.
 41 Max. load allowed for chamber interior total shelves (31 and chamber bottom, excluding load on casters mounted on product.
- casters mounted on product.

 5) Nominal value Background noise 20 dB[A].

 6) Installation of small inner door kit may effect usable
- 7) Requires sensor cover MTR-DU700SF-PW. 8) Only for MTR-5000 (data acquisition system) users. 9) MDF-DU502VH, MDF-DU502VHL, MDF-DU702VH, MDF-DU702VHL are for laboratory use.
- Appearance and specifications are subject to change

Caution: PHC Corporation guarantees this product under certain warranty conditions. However, please note that PHC Corporation shall not be responsible for any loss or damage to the contents of the product.



Preservation Equipment, Experimental Environment Equipment, Dispensary Equipment, Culturing Equipment and Drying & Sterilising Equipment for General Laboratory use

The management of the design, development, production and servicing of the above. PHC Corporation, Biomedical Division

1-1-1 Sakada, Oizumi-machi, Ora-gun, Gunma 370-0596, Japan







PHC Corporation Biomedical Division is certified for: Environmental

management system: IS014001

DISTRIBUTED BY:

