



MCO-173AIC-PE

IncuSafe

CO₂ Incubators

165 L



(option*)



(option**)



(page 2*)



Optimising cell culture outcomes and reproducibility

IncuSafe CO₂ Incubators provide precise control of CO₂ concentration and accurate, uniform, and highly responsive temperature control within the chamber. During cell culturing the inCu-saFe germicidal interior and SafeCell UV-LED lamp work continuously to prevent contamination.

Precise & Regulated Environment


InCu-saFe and SafeCell UV-LED both function to prevent contamination. The Direct Heat and Air Jacket System regulates the temperature whilst the Dual IR sensor controls the CO₂ level.

Time-Saving Decontamination

The high-speed decontamination system uses vaporised hydrogen peroxide and UV-LED light. It cleans the chamber of the incubator safely in less than three hours, achieving a minimal 6 log reduction of major contaminants.

Ease of Use & Maintenance

A new 7 inch color LCD touch panel allows full control even with gloved hands. Transfer of data is easy via a USB port. The easy-to-clean incubator interior features fully rounded corners and integrated shelf supports.

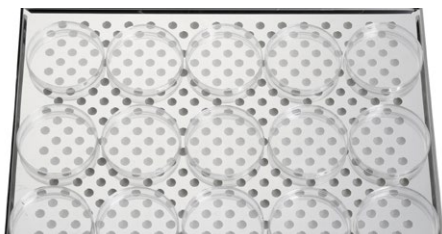

Optimum Cell Growth

Outstanding quality and performance for successful cell growth, optimal results and reproducibility. Perfect fit for the strictest and most sensitive protocols.



Efficient Workflows

Complete laboratory procedures and experiments more efficiently with less incubator downtime. Ideal for commercial applications.



Intuitive Usability

Control and visibility of the internal conditions, such as CO₂ level and temperature, is easy with the MCO-173AIC CO₂ incubator.

* Standard for Models MCO-173AICUV-PE/MCO-173AICUVH-PE
** Standard for Model MCO-173AICUVH-PE

IncuSafe CO₂ Incubators



Direct Heat and Air Jacket System

Achieves accurate, uniform, and highly responsive temperature control within the chamber, providing exceptional uniformity and rapid recovery after door-openings.

Dual IR CO₂ Sensor

The incubator's Dual IR sensor and P.I.D control enables ultra-fast CO₂ recovery without overshoot, even following multiple door-openings.


Active Background Decontamination

The exclusive inCu-saFe copper-enriched stainless steel alloy interior offers the germicidal properties of copper and the durability of stainless steel. The optional, isolated, SafeCell UV -LED lamp decontaminates circulating air and water in the humidifying pan, without harming cultured cells.

Condensation Management

The 'dew stick' -controlled by Peltier technology- condenses water on its surface, which then drips into the humidifying pan, preventing unwanted condensation in the chamber and possible contamination.

Dimensions

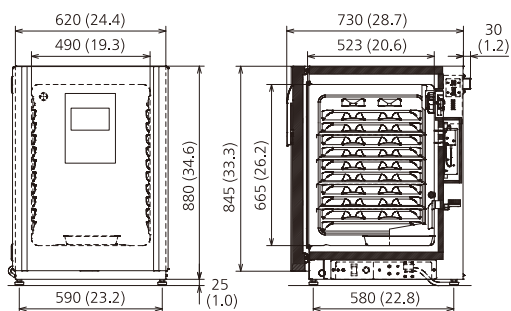

Unit: mm (inch)

#)



0123 Available in these countries as Class IIa Medical Device:
Austria, Belgium, Cyprus, France, Germany, Ireland, Italy, Liechtenstein, Luxembourg, Malta, the Netherlands, Poland, Spain, Switzerland and the United Kingdom.



Available as equipment for laboratory use in:
EEA countries, Switzerland, the United Kingdom and Turkey.

PHC Europe

A Member of PHC Group

Eikdonk 1 | 4825 AZ Breda | Netherlands
T: +31 (0) 76 543 3833

www.phchd.com/eu/biomedical

Model Number		MCO-173AIC-PE	MCO-173AICUV-PE	MCO-173AICUVH-PE
External Dimensions (W x D x H) ¹⁾	mm	620 x 730 x 905		
Internal Dimensions (W x D x H)	mm	490 x 523 x 665		
Volume	liters	165		
Net Weight	kg	69		
Performance				
Temperature Control Range & Fluctuation	°C	AT +5 ~ +50, ±0.1		
Temperature Uniformity ²⁾	°C	±0.25		
CO ₂ Control Range & Fluctuation	%	0 ~ 20, ±0.15		
Humidity Level & Fluctuation	%RH	95, ±5		
H ₂ O ₂ Decontamination		Optional	Optional	Standard
Control				
Temperature Sensor		Thermistor		
CO ₂ Sensor		Dual IR		
Display		WVGA full color LCD touchpanel		
Construction				
Exterior Material		Painted Steel (rear cover not painted)		
Interior Material		Stainless Steel Copper-Enriched Alloy		
Insulation Material		Expandable styrene-acrylonitrile beads		
Heating Method		Direct Heat & Air Jacket System		
Outer Door	qty	1		
Electric Door lock with Password		Optional	Optional	Standard
Field Reversible Door		Included		
Inner Door	qty	1 gas tight - made of tempered glass		
Shelves	qty	4 x Stainless Steel Copper-enriched Alloy		
Shelf Dimensions (W x D x H)	mm	470 x 450 x 12		
Max. Load per Shelf	kg	7		
Max. Shelf Capacity	qty	10		
Access Port	qty	1		
Access Port Position		Rear Upper Left		
Access Port Diameter	Ø mm	30		
Alarms (R = Remote Alarm, V = Visual Alarm, B = Buzzer Alarm)				
Power Failure		V-B-R		
Out of Temperature Setting		V-B-R		
High Temperature		V-B-R		
Out of CO ₂ Setting		V-B-R		
Door open		V-B		
Electrical and Noise Level				
Power Supply	V	230		
Frequency	Hz	50		
Noise Level ³⁾	dB	29		
Options				
Electrical Door Lock With Password		MCO-170EL-PW		Standard
SafeCell UV-LED® System		MCO-LUVS-PW ⁴⁾	Standard	
H ₂ O ₂ Decontamination Board		MCO-HK-PW ⁴⁾		Standard
H ₂ O ₂ Vapor Generator		MCO-50HP-PW ⁴⁾		
H ₂ O ₂ Reagent, pack of 6 bottles		MCO-20H2O2-PE		
CO ₂ Gas Pressure Regulator		MCO-010R-PW		
Automatic CO ₂ Cylinder Changeover System		MCO-21GCP-PW		
Semi-automatic one point Gas Calibration Kit		MCO-SG-PW		
InCu-saFe® Shelf		MCO-170ST-PW		
InCu-saFe® Half Tray System		MCO-25ST-PW		
Double Stacking Bracket*		MCO-170PS-PW		
Stacking Plate*		MCO-170SB-PW		
Roller Base		MCO-170RB-PW		
Communication systems ⁵⁾				
Analogue interface (4-20mA)		Standard		
Mod-bus interface board		MTR-MODB1-PW		
OPC-UA interface board		MTR-OPC1-PW		

Appearance and specifications are subject to change without notice.

¹⁾ Exterior dimensions of main cabinet only, excluding handle and other external projections.

²⁾ ±0.25°C; ambient temp 23°C - 25°C, setting 37°C, CO₂ 5%, no load

³⁾ Nominal value

⁴⁾ MCO-173AIC-PE requires MCO-HK-PW, MCO-50HP-PW and SafeCell UV-LED option for H₂O₂ decontamination

⁵⁾ Can only be fitted with one communications interface.

* If stacking two incubators, make sure the double-stacking dedicated securing hardware and spacer are used.